

# Sterisart<sup>®</sup> Family

Sterility Testing Made  
Easy and Reliable

## Benefits

- Compact, ergonomic design to cover all standard sterility testing applications
- Specially engineered needles for safe and easy piercing of stoppers and containers
- Anti-foaming inlet
- Raised sterile venting
- Tethered, large, grip-optimized filter caps
- Septum port that guarantees aseptic sampling and supplementation
- Recess in packaging that houses directly accessible outlet plugs

## Product Information

The development and manufacturing of injectable, sterile medical products under GMP conditions is one of the most challenging and sensitive issues in the (bio-)pharmaceutical industry. Before a batch of the parenterals concerned can be released by the quality control department (QC), proof of sterility must be provided by performing the sterility test according to USP <71>, Ph. Eur. 2.6.1 and JP 4.06.



## International Guidelines and Requirements

According to international pharmacopeias, parenterals (such as vaccines, mAbs, intravenous infusions, antibiotics, diabetes medication) that are injected into the human or animal body are subject to sterility testing. However, other products must be tested for sterility as well, for example:

- Ophthalmics
- Immunodiagnosics (urine | blood, etc.)
- Intermediates | APIs
- Cell culture media
- Cell banks
- Virus banks
- Medical instruments (e.g., scalpels)
- Ointments
- Creams

Various analytical methods are used depending on the volume to be analyzed; e.g., membrane filtration or direct inoculation.

Generally, the sterility of each batch of a sterile drug is verified by performing the sterility test at the latest on such pharmaceuticals packaged in their final containers. Given the role of the test method in the federal laws of some countries (e.g., U.S. CFR 610.12), this qualitative test for the presence or absence of microorganisms, yeasts and fungi also has additional, legal significance.

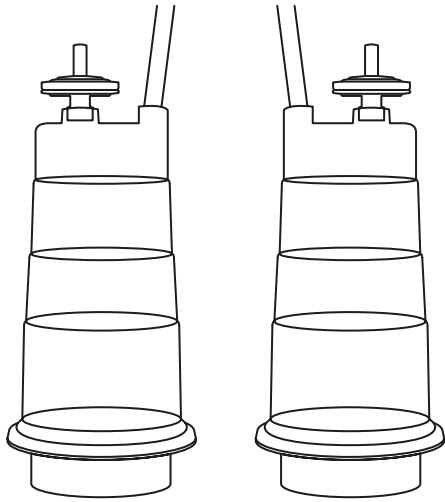
## Quality Control

As Sartorius is a recognized supplier of products and services for the pharmaceutical industry, product safety and quality are its number one priorities. To meet the quality requirements according to ISO 9001, representative samples of each lot of Sterisart® systems are subjected to destructive testing. In-process and final quality control tests further ensure high product safety.

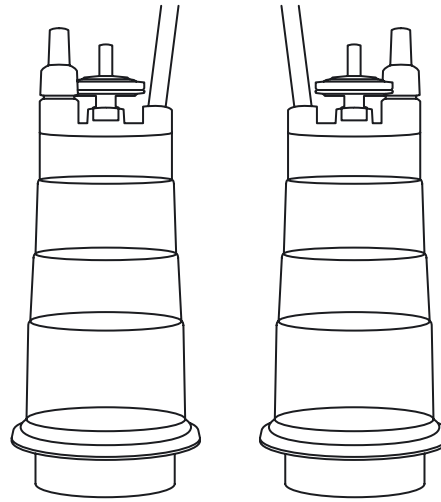
In addition to the bacteria challenge test and the Method Suitability test, the bubble point, flow rate, thickness and wetting time of the membrane are all checked.

# Technical Specifications


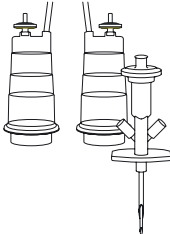
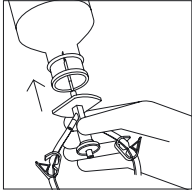

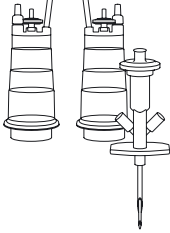
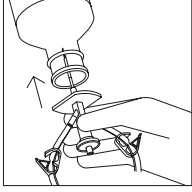
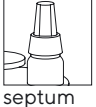

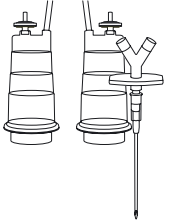
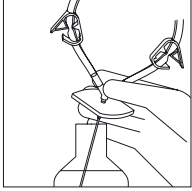
## Basic Structure of Sterisart® RC Units with a Regenerated Cellulose Membrane



Standard configuration, order no. ending in GBD

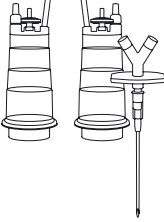
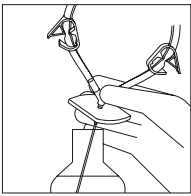

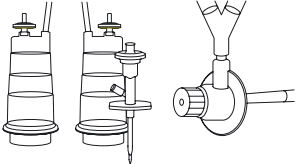
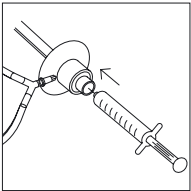

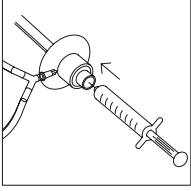

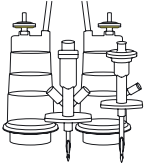
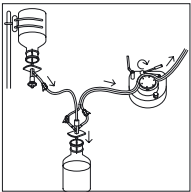
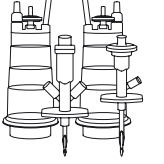
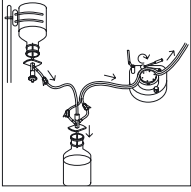

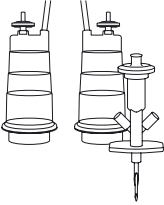
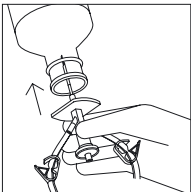
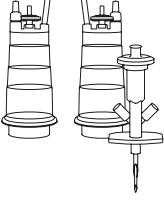
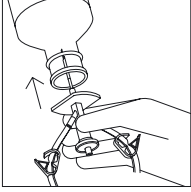



Configuration with septum, order no. ending in GSD

Sample Type	Product Container	Spike	Usage	Order Number
LVPs	Closed glass bottles with a septum 			16466-----GBD
LVPs	Closed glass bottles with a septum 		 	16466-----GSD
LVPs, SVPs	Open containers e.g. glass ampoules, glass bottles, collapsible plastic bags 			16467-----GBD

According to the United States Pharmacopeia:  
 \*LVPs: Large Volume Parenterals >100 mL  
 \*SVPs: Small Volume Parenterals <100 mL

# Basic Structure of Sterisart® RC Units with a Regenerated Cellulose Membrane

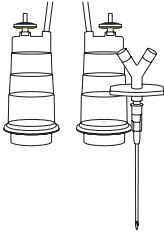
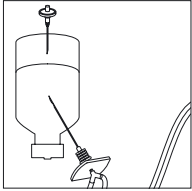
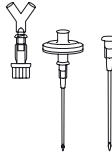
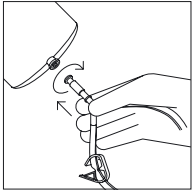
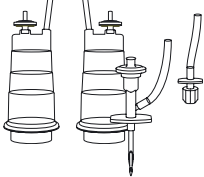
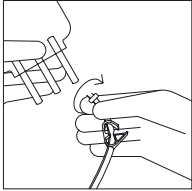
Sample Type	Product Container	Spike	Usage	Order Number
LVPs, SVPs	Open containers e.g. glass ampoules, glass bottles, col- lapsible plastic bags		  septum	16467-----GSD
SVPs	Prefilled syringes with or without a spike			16469-----GBD
SVPs	Prefilled syringes with or without a spike		  septum	16469-----GSD
Lyophilisates, soluble powders, liquid antibiotics	Closed glass bottles with a septum			16475-----GBD
Lyophilisates, soluble powders, liquid antibiotics	Closed glass bottles with a septum		  septum	16475-----GSD
SVPs	Closed glass bottles with a septum			16476-----GBD
SVPs	Closed glass bottles with a septum		  septum	16476-----GSD

According to the United States Pharmacopeia:

\*LVPs: Large Volume Parenterals >100 mL

\*SVPs: Small Volume Parenterals <100 mL

# Basic Structure of Sterisart® RC Units with a Regenerated Cellulose Membrane

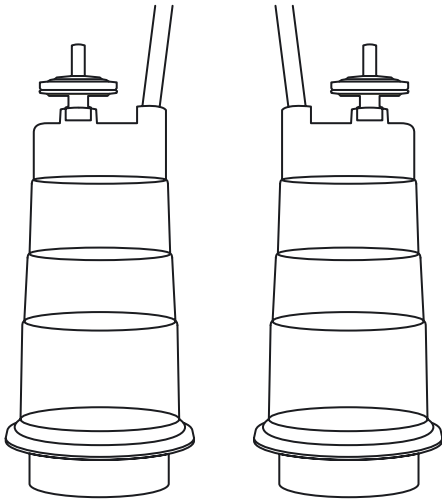
Sample Type	Product Container	Spike	Usage	Order Number
LVPs, SVPs, eye drops	Closed containers; plastic containers with blow-fill seals; e.g. bottles, ampoules			16477-----GBD
LVPs	One-connector system for testing tube assemblies and bags. Fits product containers with male Luer lock or female Luer slip connectors			16468-----GBD
LVPs	Two-connector system for testing product containers with a male Luer lock; double-needle spike for simultaneous transfer of rinsing liquid			16478-----GBD

According to the United States Pharmacopeia:

\*LVPs: Large Volume Parenterals >100 mL


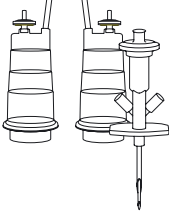
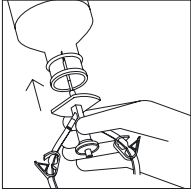

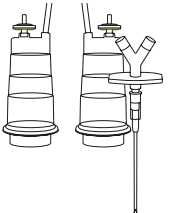
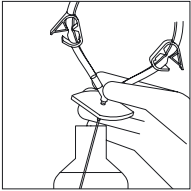
\*SVPs: Small Volume Parenterals <100 mL





Standard configuration, order no. ending in GBD

## Basic Structure of Sterisart® CA Units with a Cellulose Acetate Membrane

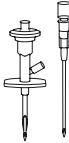

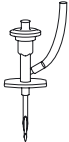

Sample Type	Product Container	Spike	Usage	Order Number
LVPs	Closed glass bottles with septum 			1646601-----GBD
LVPs, SVPs	Open containers; e.g., glass ampoules, glass bottles, collapsible plastic bags 			1646701-----GBD

According to the United States Pharmacopeia:

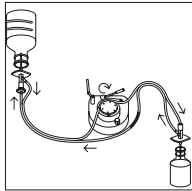
\*LVPs: Large Volume Parenterals > 100 mL

\*SVPs: Small Volume Parenterals < 100 mL

## Products for Direct Aseptic Transfer of Liquid Samples to Each Required Culture Medium

Sample Type	Product Container	Spike	Figure	Order Number
Liquids difficult to filter; medical materials	Closed glass bottles with a septum			16472-----GBD
Liquids difficult to filter; medical materials	Closed glass bottles with a septum   open glass bottles			16471-----GBD

## Sample Preparation

Sample Type	Product Container	Usage	Order Number
Powder with low solubility in closed glass bottles with a septum	Tube assembly with metal double needles of two different lengths		16470-----GBD

## Optional Accessories

Sample Type	Product Container	Usage	Order Number
Sterile venting of containers filled with rinsing solutions and culture media	Spike with 0.2 $\mu\text{m}$ sterilizing-grade filter, 4 cm, stainless steel, individually sterile-packaged, gamma-sterilized		16596-----HNK

# Applications

## Applications for Different Membranes and Systems

**Sterisart® RC** with a regenerated cellulose membrane is particularly suitable for testing aggressive, aqueous products and antibiotics.

**Sterisart® CA** with a cellulose acetate membrane has been specially designed to analyze the sterility of difficult-to-filter, viscous substances, such as emulsions, using membrane filtration.

**Transfer sets:** If the membrane filtration method cannot be used for sterility testing of liquid products, the Sartorius sterile transfer sets can be used for aseptic transfer of these products to liquid culture media for performing the direct inoculation method.

## Overview of the Sterisart® Systems

Gamma-sterilized systems with regenerated cellulose (RC) membrane  
16466-GBD, 16467-GBD, 16468-GBD, 16469-GBD, 16470-GBD, 16471-GBD, 16472-GBD, 16475-GBD, 16476-GBD, 16477-GBD, 16478-GBD,

Gamma-sterilized systems with a septum and a regenerated cellulose (RC) membrane  
16466-GSD, 16467-GSD, 16469-GSD, 16475-GSD, 16476-GSD

Gamma-sterilized systems with a cellulose acetate (CA) membrane  
1646601-GBD; 1646701-GBD

### Note:

The primary packaging of the systems as well as the primary packing of the transfer kits are gas-tight. This enables them to be used directly in isolators and prevents unnecessary rinsing steps.

### EXPAND® Training

Available on request in your company or in our Training center.





# Sterisart® Universal Pumps for Transfer of Liquids to the Sterisart® Units

**16420 with a display and an integrated barcode scanner**  
**16419 Basic version**

Order number	Description
1Z---0004	Sterisart® Easy Configuration Software Drag & drop software, designed to increase process reliability for sterility testing
1ZG---0023	Drainage container cover for Sterisart® sterility test systems
1ZE---0033	Trolley for Sterisart® Universal Pump
1ZGF---0020	Tray for 10 Sterisart® units
1ZE---0039	Trolley for pump
1ZA---0002	Drain tubing
1ZG---0028	Drainage container and cover for Sterisart® Sterility units
1ZGL---0033	Bottle holder without wing nut
1ZF---0007	Wing nut

Order number	Description
1ZGD---0031	Stainless steel cover for the rotor
1EE---0010	External barcode scanner
1ZE---0050	Isolator installation kit
1ZG---0024	Drainage container cover for competitor's sterility test consumables
1ZG---0014	Adapter for Sterisart® systems in Equinox pumps; pkg. of 2
1ZA---0028	Drainage container for Sterisart® pump 16419   16420

## Service

EXTEND® instrument services for Sterisart® Universal pumps

- Installation qualification and operational qualification (IQ | OQ)
- Preventive maintenance



# General Technical System Details and Regulatory Requirements

The Sterisart® systems comply with GMP requirements with respect to sterility testing as well as to Regulation (EC) No. 1907/2006 (REACH) with particular reference to the plasticizer DEHP; ED/108/2014, ED/ 67/2008.

## Sterilization

Gamma-sterilized at 25 kGy in compliance with DIN EN 552 and ISO 11137

Sterilization indicator on each carton; color change: from orange to red

## Shelf Life

Gamma-sterilized units:

Sterile for up to 3 years after the date of manufacture;  
guaranteed for a minimum of 6 months after delivery



## Dimensions and Weight

Box (W×D×H) 28.1×27.3×24.6 cm

Weight of carton with 10 Sterisart® units Approx. 2.2 kg

Gas-tight individual packaging (W×D×H) 5.5×13×27 cm

Weight of Sterisart® unit with packaging Approx. 185 g

From bottom to sterile air filter red cap attached (H×D) Approx. 13.8×5.7 cm

Weight of the Sterisart® unit without primary packaging Approx. 150 g, depending on the system version

## Packaging

Primary packaging material OPA | PE sealing foil

Transparent film A-PET/PE

Transport packaging material Polyamide (PA) and polyethylene (PE)

Transparent film PETG (polyethylene terephthalate)



## Sterisart® System

Method Suitability tests are performed on Sterisart® systems in compliance with the international pharmacopeias; USP <71> and Ph. Eur. 2.6.1.; an extractables profile has been created in compliance with the USP 23 and Ph. Eur. 3. The detection limits are below the requirements as specified for “water for injection”.

Equal distribution of the product to the Sterisart® containers with a max. validated deviation of 10 %

Regenerated cellulose (RC) membrane	
Pore size	0.45 µm nominal; specified according to international pharmacopeias; USP <71>; Ph. Eur. 2.6.1
Effective filter area	15.7 cm <sup>2</sup>
Integrity test	2.5 bar
Membrane thickness	Approx. 150 – 170 µm
Transparent film	PETG (polyethylene terephthalate)

Cellulose acetate (CA) membrane	
Pore size	0.45 µm nominal; specified according to international pharmacopeias; USP <71>; Ph. Eur. 2.6.1
Effective filter area	14.5 cm <sup>2</sup>
Integrity test	1.5 bar
Membrane thickness	115 – 145 µm
Transparent film	PETG (polyethylene terephthalate)

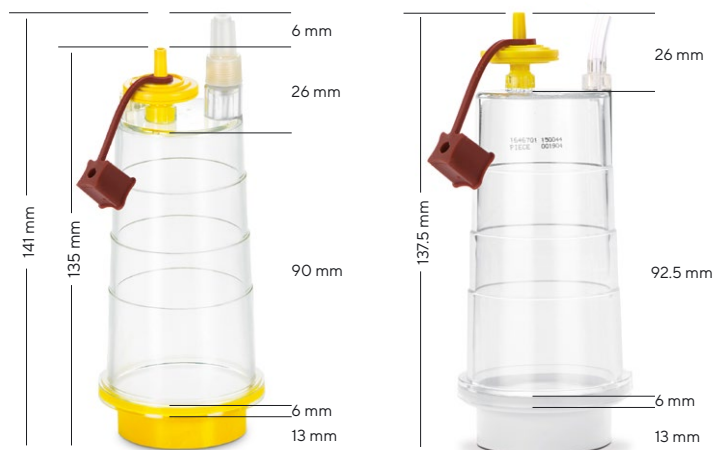
Hydrophobic (sterile venting) filters	
Membrane	0.2 µm polytetrafluoroethylene (PTFE); validated according to HIMA for the retention of <i>Brevundimonas diminuta</i>
Burst pressure of the sterile air filter	At least 6 bar
Housing	Acrylic-based multipolymer
Water permeability pressure   penetration pressure	> 3 bar
Transparent film	PETG (polyethylene terephthalate)

Sterisart® container	
Upper part   lower part	Styrene acrylonitrile (SAN)
Burst pressure of the housing	> 5 bar
Max. operating pressure	3 bar at 20 °C
Max. operating temperature	50 °C
Capacity	120 mL (50 mL, 75 mL and 100 mL graduated marks)
Red caps	Silicone
Integral membrane	Membrane incorporated in the system by a special clamping technology
Syringe holder	Styrene acrylonitrile (SAN) Sterisart® system, 16469
Female Luer lock and Luer slip connectors	Acrylonitrile butadiene styrene (ABS) Sterisart® system 16478



Accessory Materials Kit	
Wing nuts	Polyethylene (PE)
Tubing and Spikes	
Tubing material and length	PVC, 80 cm; Additionally in available as silicone tubing, order no. 16469; 60 cm
Sampling needle	Polycarbonate and stainless steel (Sterisart® system 16467)
Double-needle spike	Acrylonitrile butadiene styrene (ABS) and stainless steel
Needle for sterile venting	Polypropylene (PP) and stainless steel; PTFE membrane; Housing made of methyl acrylate-butadiene-styrene (MBS)

	Length	Outer Diameter
Needle for sterile venting of Sterisart® systems 16467, 16477 and 16468	40 mm	1.6 mm
Needle for Sterisart® system 16467	52 mm	1.5 mm
Needle for Sterisart® system 16468	60 mm	1.5 mm
Double-needle spike (long) for Sterisart® system 16466	41 mm	2.8 mm
Double-needle spike (short) for Sterisart® system 16476	21 mm	2.8 mm
Double-needle spike for Sterisart® systems 16469, 16471, 16475, 16478	35 mm	2.8 mm
Double-needle spike for Sterisart® system 16475	23 mm	2.8 mm
Double-needle spike (short) for Sterisart® system (one connector) 16472	22 mm	2.8 mm
Double-needle spike (short) for Sterisart® system (one connector) 16470	23 mm	2.8 mm
Double-needle spike (long) for Sterisart® system (one connector) 16470	37 mm	2.8 mm



#### Septum Material (only in systems ending in GSD)

Protective cap	Polyethylene
Septum material	Polysoprene and acrylonitrile butadiene styrene (ABS)

**000 «Диаэм»**

**Москва**  
ул. Магаданская, д. 7, к. 3 ■ тел./факс: (495) 745-0508 ■ sales@dia-m.ru

**www.dia-m.ru**

**С.-Петербург**  
+7 (812) 372-6040  
spb@dia-m.ru

**Новосибирск**  
+7(383) 328-0048  
nsk@dia-m.ru

**Воронеж**  
+7 (473) 232-4412  
vrn@dia-m.ru

**Йошкар-Ола**  
+7 (927) 880-3676  
nba@dia-m.ru

**Красноярск**  
+7(923) 303-0152  
krsk@dia-m.ru

**Казань**  
+7(843) 210-2080  
kazan@dia-m.ru

**Ростов-на-Дону**  
+7 (863) 303-5500  
rnd@dia-m.ru

**Екатеринбург**  
+7 (912) 658-7606  
ekb@dia-m.ru

**Кемерово**  
+7 (923) 158-6753  
kemerovo@dia-m.ru

**Армения**  
+7 (094) 01-0173  
armenia@dia-m.ru

